

Freezers - Ultrafreezers Platilab & Irilab Series





- → **PLATILAB**: top security and performances with reliable temperature control.
- → IRILAB: combining the high reliability of PLATILAB with the possibility to store up to 60.000 vials in less than one square meter thanks to V.I.P. insulation.

-85.0°C



Platilab - Irilab Series /Optional





→ Chest and up-right freezers at -40°C / -86°C for freezing and long term storage of biological samples.

- \rightarrow Low energy consumption.
- \rightarrow V.I.P. Technology for footprint optimization (Irilab models).
- → Microprocessor control board with seven independent sections for top reliability in case of failures (NEXT models).
- \rightarrow Alarm logs in text and graphs.
- \rightarrow Data export on SD memory card.
- \rightarrow Access control.
- → Double door models with two indipendent chambers (NEXT DD models).





 \rightarrow Professional line.

- \rightarrow Models **H** and **V** Working temperature: -86°C.
- \rightarrow Models SH and SV Working temperature: -40°C.

Technical features

Structure /Fully metallic in heavy-gauge phosphate-coated, powder-painted steel plate with interior in top quality stainless steel AISI 304.

Both up-right and chest models are installed on castors and adjustable feet. **Door** /PLATILAB **STD** models equipped with mechanical lock and key;

IRILAB **STD** models equipped with magnetic lock and key.

In **PLUS** version the door is equipped with patented transponder system **BIOGUARD**[®], which enables the door opening by a card to authorized personnel only.

The system guarantees a complete access control. In case of power outage, the buffer battery maintains the system on for 24 hours.

Heated pressure compensation valve allows an easy and quick opening of the door. **Port hole** /Diameter 23 mm, with rubber cap for the insertion of additional probes. **Inner doors (only for up-right models)** /

PLATILAB models: antishock polystyrene panels with magnetic lock and special gaskets. IRILAB models: composite materials (aluminum - pvc - aluminum) panels and magnetic stop. Shelves (only for up-right models) / Stainless steel adjustable in height. Insulation /

PLATILAB models: thickness 140 mm. High pressure (40 kg/mc) polyurethane foam. IRILAB models: thickness 70 mm. Vacuum Insulated Panels (V.I.P.) and polyurethane foam. **Cooling system** /Hermetically sealed, two stages (Cascade cycle) compressor (-86°C) and single stage compressor (-40°C). Refrigerant HCFC-free.

Adjustment and control system /PLATILAB e IRILAB models: microprocessor controller E2003, with two independent sections, the first one to set and control temperature, the second one to manage the alarm system and display temperature.

Rechargeable buffer battery (12 V, 2Ah) to maintain the alarm system on for 48 hours in case of power outage. Configured for a remote replication of alarms.

01 02 03









Next models add to Platilab - Irilab freezers enhanced display and data logging features by means of **Cold Brain™**, the superior microprocessor controller with seven independent sections.

Microprocessor controller **Cold Brain™**, with seven independent sections to maintain the proper working conditions even in presence of failures.

The multiple high quality sensors (four PT100 and four NTC) and the efficient and flexible alarm system guarantee the maximum security in sample control.

Cold Brain[™] key features

- **Multi language** compliant to European rules.
- Wide graphic display (240x128pixel), with regulation and alarm sections.
- **Touch keyboard** using 8 soft keys with context dependent functions.
- On-Screen System Messaging: displays set-points, details of alarms, door openings and operating conditions.
- **Integrated Data Recorder:** traditional paper chart recorder is replaced by electronic recorder. Data (alarms and parameters) are stored for 3 months at a sample interval of 1 min.
- Acoustic and Visual Alarms for high and low temperature conditions, power failure, low battery, malfunctions and filter cleaning.
- Alarm Log provided via text and graphis, to record door openings, power failure, electronic fault, internal probe and ambient temperature probe failures, fan and condenser malfunctions.
- **Back up battery** (up to 48 hours): maintains settings and activates visual and acoustic alarms during power outage.
- Battery charge indicator.
- **Prearranged for CO2 & LN2 Backup Systems:** simplifies and makes cheaper the installation.
- Remote Alarm Contacts provides connection to centralized monitoring systems.
- Secure Digital (SD) card to download data on PC.
- Ethernet Interface (Optional Wireless) to connect to a laboratory pC network. Through Cryolog2 software, it is possible to monitor parameters and alarms of multiple freezers (up to 30).
- **Protected On-Off Switch** to prevent accidental shut off.
- Low noise.

04 05



01. E2003 with Bioguard[®].

- 02. Irilab 800 with 4 inner doors.
- 03. Platilab 500 with 3 inner doors and racks.
- 04. Irilab NEXT 800.
- **05.** COLD BRAIN[™] detail.

Double door Freezers: Next™ DD Models

Indipendent chambers /Two separated chambers, accessible through independent doors. **Structure** /Heavy-gauge phosphate-coated, powder-painted steel plate with interior in top quality stainless steel AISI 304. Installed on castors and adjustable feet, for easy movement. All our freezers fit with standard doors.

Port hole /Diameter 23 mm, with rubber cap for the insertion of additional probes. **Door /STD** models: equipped with magnetic lock and key.

The **PLUS** version door is equipped with patented transponder system **BIOGUARD**[®], which enables the door opening by a card to authorized personnel only. The system guarantees a complete access control.

In case of power outage, the buffer battery maintains the system on for 24 hours. Heated pressure compensation valve allows an easy and quick opening of the door. **Inner doors** /Panels in composite material (aluminum–pvc–aluminum) and magnetic stop.

Shelves (only for up-right models) /Stainless steel adjustable height. Insulation /Thickness 70 mm. Vacuum Insulated Panels (V.I.P.) and polyurethane foam. Cooling system /Double stage hermetic system. HCFC-free refrigerant.Adjustment and control system /Microprocessor controller Cold Brain[™], with seven independent sections to maintain the proper working conditions even in presence of failures. The multiple high quality sensors (four PT100 and four 4 NTC) the efficient and flexible alarm system quarantee the maximum security in sample control.

Cold Brain[™] key features

- **Multi language** compliant to European rules.
- Wide graphic display (240 x128 pixel), with regulation and alarm sections.
- **Touch keyboard** using 8 soft keys with context dependent functions.
- **On-Screen System Messaging**: displays set-points, details of alarms, door openings and operating conditions.
- Integrated Data Recorder: traditional paper chart recorder is replaced by electronic recorder. Data (alarms and parameters) are stored for 3 months at a sample interval of 1 min.
- Acoustic and Visual Alarms for high and low temperature conditions, power failure, low battery, malfunctions and filter cleaning.
- **Alarm Log** in text and graphs, to check door openings, power failure, electronic fault, internal probe and ambient temperature probe failures, fan and condenser malfunctions.
- **Backup battery** (up to 48 hours): maintains settings and activates visual and acoustic alarms during power outage.
- Battery charge indicator
- Prearranged for CO2 & LN2 Backup Systems: simplifies and makes cheaper the installation.
- **Remote Alarm Contacts** provide connection to centralized monitoring systems.
- Secure Digital (SD) card to download data on PC.
- Ethernet interface (Optional wireless) to connect to a laboratory PC network. Through Cryolog2 software, it is possible to monitor parameters and alarms of multiple freezers (up to 30).
- Protected On-Off Switch to prevent accidental shut off.
- Low noise.

06 07





06. Irilab NEXT 520 DD.07. Irilab NEXT 520 DD with doors opened.

Technical specifications

	Horizontal Models	Horizontal Models						
Model	PLATILAB 110		PLATILAB 370		PLATILAB 550			
	SH H		SH H		SH H			
Nominal capacity (I)	110		3	70		55	i0	
Temperature range (°C)	-20 /-45	40 / -86	-20 / -45	-40 / -86	-20	/ -45	-40 / -86	
Supply voltage (V)*		230 V /50 Hz - 115 V /60 Hz - 230 V /60 Hz						
Refrigerant	R404 mixe	R404 mixed refrigerant		R404 mixed refrigerant		R404 mixed refrigerant		
External dimensions (mm)	730 × 850 × 108	30	1780 × 855 × 1140		2350 × 855 × 1140		i5 × 1140	
nternal dimensions (mm)	450 × 560 × 45	450 × 560 × 450		1050 × 520 × 670		1590 × 520 × 670		
Max Boxes capacity (n°)	63	63		210		330		
Overall weight with standard packaging (Kg)	175	205	274	299	3	24	360	
Noise level (dbA)	52							
	Vertical Models							
Model	PLATILAB 340		PLATILAB 500		IRILAB 800			
	SV V		SV V		SV V			
Nominal capacity (I)	340	D 5		00		800		
Femperature range (°C)	-20 /-45	40 / -86	-20 / -45	-40 / -86	-20	/ -45	-40 / -86	
Supply voltage (V)*			230 V /50 Hz - 115 V	/60 Hz - 230 V /60 Hz	Z			
Refrigerant	R404 mixe	d refrigerant	R404	mixed refrigerant	R404 mixed refrigerant			
External dimensions (mm)	765 × 900 × 198	35	1015 × 90	00 × 1985	1086 × 890 × 1995			
nternal dimensions (mm)	450 × 530 × 135	450 × 530 × 1350		700 × 530 × 1350		906 × 609 × 1472		
Max Boxes capacity (n°)	255		420		600			
Overall weight with standard packaging (Kg)	243	260	293	330	3	50	390	
Noise level (dbA)			<52				<53	
	Horizontal Models Vertical Models							
Model	PLATILAB Next 370	PLATILAB Next 370 PLATILAB Next 5		PLATILAB Next 340		PLATILAB Next 500		
Nominal capacity (I)	370	370		550 340		500		
Femperature range (°C)		340						
upply voltage (V)*		230 V /50 Hz - 115 V /60 Hz - 230 V /60 Hz						
Refrigerant	mixed refrigerant							
external dimensions (mm)	1780 × 855 × 1140	2350	0 × 855 × 1140	765 × 900 × 1985		1015	× 900 × 1985	
nternal dimensions (mm)	1050 × 520 × 670	1050 × 520 × 670 1590		450 × 530 × 1	1350 700 × 530 × 1350		× 530 × 1350	
Max Boxes capacity (n°)	210	210		330 252		420		
Overall weight with standard packaging (Kg)	299	360		260			330	
Noise level (dbA)			5	2				
	Vertical Models							
Model	IRILAB Next 520	IRILAB Next 520			IRILAB Next 800			
	V	V - DD		V		V - DD		
Nominal capacity (I)	· · · · · · · · · · · · · · · · · · ·	520				800		
Femperature range (°C)			-40	/-86				
Supply voltage (V)*		230 V /50 Hz - 115 V /60 Hz - 230 V /60 Hz						
Refrigerant				frigerant				
External dimensions (mm)	7305	730×854×1980 1086×890×1995						
nternal dimensions (mm)		586×613×1470 906×609×1472						
Max Boxes capacity (n°)	384	1			600 / 432			
nan bones capacity (in)	1 504		200					
Overall weight with standard packaging (Kg)	260		274	390			407	

 $\stackrel{>}{\rightarrow}$

Illustration, description and technical data may be changed without notice. -40°C and -86°C freezers certified according to 93/42/ECC directive are also available (please refer to **Blood Bank Refrigerators and Freezers** catalogue). (*) 230V/60Hz and 115V/60Hz models are also available.





Angelantoni Life Science S.r.l.

Località Cimacolle, 464 - 06056 Massa Martana (PG) Italy ph. (+39) 075.89551 (a.r.) - fax (+39) 075.8955312 biomedical@angelantoni.it

www.angelantonilifescience.it

